# KÖNIG PHOTO

## KN-MICRO700

**MANUAL (p. 2)  
STEREO MICROSCOPE**

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**MANUALE (p. 10)  
MICROSCOPIO STEREO**

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## ENGLISH

### NOTICE

1. A microscope is a precision instrument, so handle with care, avoiding impact or abrupt movement during transportation.
2. Microscope ought to be placed in a dry and clean place. Do not expose the microscope directly to the sun. Avoid high temperature and violent vibration. Following environment is required: Indoor temperature: 0° - 40°, Max relative humidity: 85%.
3. To keep the image clear, do not leave fingerprints or stains on the surfaces of the lens.

### Maintainance

1. All lenses must be kept clean. Fine dust on the surface of the lens should be blown off with hand blower or wiped off gently with a soft lens tissue. Fingerprints or oil marked on it should be wiped off with a tissue moistened with small amount of 3:7 mixture of alcohol and ether.
2. Never use the organic solution to clean the other surface (especially the plastic surfaces). If necessary, please choose the neutral detergent.
3. Do not take the microscope apart if you think that it is damaged.
4. After using, cover the microscope with the dust-cover provided and store it in a dry and clean place free from moisture to prevent rust.

### Replace the lamp

1. Before changing the lamp, you must cut off the power and wait until the lamp has cooled down.
2. When changing the top lamp, loosen the lock screw on the lamp house and take down the lamp cover and replace it with a new one.
3. When changing the bottom lamp, remove the stage from the base and then replace it with a new one.

### Power supply:

220 V-240 V $\pm$ 10%, 50-60 Hz

Lamp: Tungsten bulb 12 V/10 W or Halogen bulb

12 V/10 W

**Declaration of conformity**

We,  
Nedis B.V.  
De Tweeling 28  
5215MC 's-Hertogenbosch  
The Netherlands  
Tel.: 0031 73 599 1055  
Email: info@nedis.com

Declare that product:

Name: KONIG  
Model: KN-MICRO700  
Description: Microscope

Is in conformity with the following standards

EMC: EN61010-1/A2/07.95  
LVD: EN61010-1/04.93

Following the provisions of the 73/23/EEC directive.

's-Hertogenbosch, 2007, December 12

Mrs. J. Gilad  
Purchase Director

**Safety precautions:**

To reduce risk of electric shock, this product should ONLY be opened by an authorized technician when service is required. Disconnect the product from mains and other equipment if a problem should occur. Do not expose the product to water or moisture.

**Maintenance:**

Clean only with a dry cloth. Do not use cleaning solvents or abrasives.

**Warranty:**

No guarantee or liability can be accepted for any changes and modifications of the product or damage caused due to incorrect use of this product.

**General:**

Designs and specifications are subject to change without notice.

All logos brands and product names are trademarks or registered trademarks of their respective holders and are hereby recognized as such.

**Attention:**

This product is marked with this symbol. It means that used electrical and electronic products should not be mixed with general household waste. There is a separate collections system for these products.

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Copyright ©